

Plasmid Midi Kit

Product #: Bi2M-PlasMidi

Quantity: 10 reactions

Product Description

- Designed for rapid and cost-effectives medium-scale preparation of high quality plasmid DNA from recombinant *E.coli* cultures.
- Based on modified SDS-alkali lysis method and utilizes an exclusive silica-based membrane technology in the form of a convenient spin column. The kit utilizes proprietary glass fiber membrane for specific plasmid adsorption and a specialized buffer combination.
- Effectively extracts various endotoxin-free plasmids (≤ 0.1 EU/µg endotoxin content) from cultured bacteria.
- $\circ~$ Typical plasmid yield ranges from 100 to 400 $\mu g.$
 - The actual plasmid yield will depend on the plasmid copy number & culture density.
- The purified plasmid DNA can be used in all molecular biology procedures such as conventional digestion with restriction enzymes, PCR, transformation and sequencing.

Storage

- Store RNaseA at -20 °C. All other items can be stored at room temperature for up to 12 months. For longer storage, store at 4°C.
- After addition of RNase A, the resuspension solution **Buffer P1** should be stored at 4°C.

Kit Components	
Buffer BL	25 mL
Buffer P1	30 mL (<i>note: add 600 μL RNase A, and store at 4 °C</i>)
Buffer P2	30 mL (note: seal tightly after each use)
Buffer P3	30 mL (<i>note: pre-cool to 4 ^oC before use</i>)
Buffer ER	10 mL
Buffer ED	40 mL
Buffer PW	16 mL (note: add 64 mL 100% ethanol per bottle before use)
Buffer TE	12 mL (note: pre-heat to 60 °C before use)
RNase A	600 μL

Column Filters	10 columns
HiBind DNA Midi Column	10 columns
Collection Tube	10 tubes

Important Preparatory Notes (PLEASE READ CAREFULLY)

- 1. Prepare a 42°C water bath or heating module.
- 2. Add the entire **RNase A** provided to **Buffer P1** before use; it can then be stored at 4°C for up to 6 months.
- 3. Add 64 mL of anhydrous ethanol to Buffer PW before use.
- 4. If **Buffer P2** has precipitated, heat it in a 37°C water bath for a few minutes until it becomes clear. Close the lid immediately after use to prevent prolonged exposure to air.
- 5. Pre-cool Buffer P3 at 4°C before use.
- 6. Always keep buffer bottles tightly closed and keep columns sealed in the enclosed zip-lock bag.
- 7. Proper laboratory safety practices should be employed, including the use of lab coats, gloves, and eye protection.

Protocol

- 1. **Column Preparation:** Add 2 mL of **Buffer BL** to the **HiBind DNA Midi Column**, centrifuge at **4,500×g** for **2 minutes** at **room temperature**. Discard the filtrate and return the HiBind DNA Midi Column to the Collection Tube. It is recommended to use the prepared column immediately.
- Bacterial Culture Harvest: Harvest bacterial culture from 20-50 mL of fresh bacterial culture incubated for 12-16 h by centrifugation at 8,000 x g for 5 min at room temperature. Decant the supernatant and remove all remaining medium.

- Note: For **high-copy-number plasmids**, do not process more than 50 mL of bacterial culture. If more than 50 mL, the Hi-Bind DNA Mini Column capacity will be exceeded and no increase in plasmid yield will be obtained.
- Note: For **low-copy-number plasmids**, it may be necessary to process larger volumes of bacterial culture (up to 100 mL) to recover sufficient quantity of DNA.
- 3. **Resuspend Bacteria:** Add **2.5 mL** of **Buffer P1** (*note: ensure RNase A is added*) and thoroughly resuspend the bacterial pellet using a pipette or vortex. Ensure the bacteria are fully dispersed to optimize lysis and achieve high-quality plasmid extraction.
- 4. **Cell Lysis:** Add **2.5 mL** of **Buffer P2**, then gently invert the tube 8–10 times to lyse the cells completely. <u>Avoid vigorous shaking</u> and complete this step within **5 minutes**. The solution should become clear and viscous. If not, the culture density may be too high; consider reducing the amount of bacteria.
- Neutralization: Add 2.5 mL of pre-chilled Buffer P3, then gently invert the tube 10–12 times until a precipitate forms. Centrifuge at 8,000×g for 15 minutes at room temperature. Carefully pour the supernatant into the Column Filter, filter by pushing the handle, and collect the filtrate in a nuclease-free 15 mL centrifuge tube (not provided).
- Endotoxin Removal: Add 750 μL of Buffer ER and mix by inverting 3–5 times.
 Place the solution on ice for 10 minutes, during which the solution should turn transparent blue.
- 7. Incubate and Separate Layers: Incubate at 42°C for 5 minutes until the solution becomes turbid. Centrifuge at 4,500×g for 5 minutes at room temperature. The solution will separate into a clear upper aqueous layer and a blue lower layer. Carefully transfer the upper layer to a new nuclease-free 15 mL centrifuge tube.
- 8. Ethanol Precipitation: Add anhydrous ethanol equal to half the volume of the collected solution. Mix by inverting and transfer to the HiBind DNA Midi Column (maximum 4 mL).

EXAMPLE 1 IN THE STATE OF THE

- Column Washing: Centrifuge at 4,500×g for 3 minutes at room temperature. Discard the filtrate and return the HiBind DNA Midi Column to the Collection Tube. Repeat the loading step if necessary.
- 10. First Wash with Buffer ED: Add 3.5 mL of Buffer ED to the HiBind DNA Midi Column and centrifuge at 4,500×g for 3 minutes at room temperature. Discard the filtrate and return the column to the Collection Tube.
- First Wash with Buffer PW: Add 3.5 mL of Buffer PW (*note: ensure anhydrous ethanol has been added*) to the HiBind DNA Midi Column, then centrifuge at 4,500×g for 3 minutes at room temperature. Discard the filtrate and return the column to the Collection Tube.
- 12. Second Wash with Buffer PW: Repeat Step 11.
- Final Spin: Centrifuge at 4,500×g for 10 minutes at room temperature. Place the HiBind DNA Midi Column into a <u>new nuclease-free 15 mL centrifuge tube</u>. Open the lid and let it sit for 10 minutes at room temperature to allow complete ethanol evaporation.
- Elution: Add 100–500 μL of Buffer TE or endotoxin-free water to the center of the column membrane and let it sit at room temperature for 5 minutes. Centrifuge at 4,500×g for 5 minutes.
 - For higher plasmid yield, reapply the eluted solution to the HiBind DNA Column, allow it to sit for 5 minutes, and centrifuge again at 4,500×g for 5 minutes.
- 15. **Storage:** Store the purified plasmid DNA at -20°C or or -80°C for long-term preservation. Avoid repeated freeze-thaw cycles.

NOTES:

- 1. DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, MEDIRES CORP. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.
- 2. Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of this product, you accept the terms and conditions of all applicable limited Use Label Licenses.
- 3. ©2024 MEDIRES CORP. All rights reserved. All trademarks are the property of MEDIRES CORP. and its subsidiaries.
- 4. For Research Use Only. Not for use in diagnostic procedures.