

## Genomic DNA Extraction Kit (Spin Column Based)

**Product #:** G3633-50T

**Quantity:** 50 reactions

### Product Description

- Designed for rapid and cost-effective purification of genomic DNA from animal tissues, mammalian cells, microbial cells and whole blood.
- High purity genomic DNA extraction is achieved by combining a specially optimized lysis buffer with a proprietary silica-based membrane technology in the form of a convenient spin column.
- Rapidly extracts and purifies 10-30 µg of high purity genomic DNA from:
  - 2-30 mg of animal tissue
  - $10^6$  to  $10^7$  freshly cultured cells
  - 5-100 µL of whole blood (including anticoagulant) with non-nucleated or nucleated red cells
- Genomic DNA obtained is ready for applications such as PCR, Southern blotting, RAPD, AFLP, RFLP, sequencing and other molecular experiments.

### Kit Components – BAG #1 (store at room temperature)

Buffer GA	12 mL
Buffer GA2	25 mL x 2
Buffer GB	12 mL
Buffer PD	12 mL ( <i>note: add 18 mL 100% ethanol per bottle before use</i> )
Buffer PW	24 mL ( <i>note: add 56 mL 100% ethanol per bottle before use</i> )
Buffer TE	10 mL
DNA Spin Columns	50 columns
Collection Tubes	50 columns

### Kit Components – BAG #2 (store at -20°C)

Proteinase K	1 mL
RNase A	200 µL

### Additional Required Reagents

100% Ethanol

**Important Preparatory Notes (PLEASE READ CAREFULLY)**

1. If precipitate forms in **Buffer GA**, **Buffer GA2** or **Buffer PD** during storage, solution should be re-dissolved by incubating at 37°C, then cooled back down to 25°C before use.
2. Add **18 mL 100% ethanol** to the **Buffer PD** and mark the bottle on the label.
3. Add **56 mL 100% ethanol** to the **Buffer PW** and mark the bottle on the label.
4. Always keep buffer bottles tightly closed and keep columns sealed in the enclosed zip-lock bag.
5. Proper laboratory safety practices should be employed, including the use of lab coats, gloves, and eye protection.

**RNA Extraction Protocol****1. Sample preparation:**

- **Animal tissue:** Transfer 2-30 mg of fresh or cryopreserved animal tissue to an appropriately sized Nuclease-free tube. Add two or three 4 mm stainless steel grinding beads (Cat# G0104) per sample. Immediately add **200 µL** of **Buffer GA**. Use an automated homogenizer (TissueLyser) to homogenize thoroughly (1 to 3 min @ 20-50 Hz). Let the sample stand at room temperature for 2 minutes.
- **Adherent cells:** Remove the growth medium from the cells and wash cells with 1×PBS buffer (pH 7.4). Harvest  $10^6$  to  $10^7$  cells and transfer to an appropriately sized Nuclease-free tube. Discard the supernatant. Immediately add **200 µL** of **Buffer GA** and mix well by vortex. Let the sample stand at room temperature for 2 minutes.
- **Non-adherent suspension cells:** Transfer  $10^6$  to  $10^7$  cells to an appropriately sized Nuclease-free tube and centrifuge at 1,000 x g for 5 minutes at 4°C to pellet. Discard the growth medium from the tube. Add

**200 µL of Buffer GA** and mix well by vortex. Let the sample stand at room temperature for 2 minutes.

- **Microbial cells:** Grow a fresh 10 mL overnight culture. After incubation, centrifuge the culture at 12,000 x g for 10 minutes to harvest the cells. Discard the supernatant and resuspend the pelleted cells in **1 mL of Buffer GA2** (cell pellet can be stored at -80°C before addition of Buffer GA2 as a safe stopping point). Transfer the cell suspension to a 1.5 mL microcentrifuge tube containing 100 mg of 0.1 mm acid-washed glass grinding beads (Cat# G0310) for bacterial samples, and 100 mg of 0.5 mM acid-washed glass grinding beads (Cat# G0350) for yeast/fungal samples. Homogenize thoroughly using an automated homogenizer (TissueLyser) for 3 minutes at 30 Hz. Transfer 200 µL of the homogenized sample to a new 1.5 mL microcentrifuge tube and discard the remainder. Add **200 µL of Buffer GA** to the sample, mix well by vortexing, and let the sample stand at room temperature for 2 minutes.
- **Blood (non-nucleated erythrocytes):** Transfer 50-100 µL whole blood of non-nucleated erythrocytes (including anticoagulant) to an appropriately sized Nuclease-free tube. Add **200 µL of Buffer GA** and mix well by vortex. Let the sample stand at room temperature for 2 minutes.
- **Blood (nucleated erythrocytes):** Transfer 5-20 µL whole blood of nucleated erythrocytes (including anticoagulant) to an appropriately sized Nuclease-free tube. Add **200 µL of Buffer GA** and mix well by vortex. Let the sample stand at room temperature for 2 minutes.

## **2. Preparation and Digestion:**

- Add **20 µL of Proteinase K** and **4 µL of RNase A** to the sample from **step 1**.
- Incubate at 56°C for 30 minutes, mixing every 10 minutes to accelerate tissue lysis.
- **Note:** Extend incubation time for materials that are difficult to lyse.

**3. Centrifugation:**

- Centrifuge at 12,000 x g for 2 minutes.
- Transfer the supernatant to a nuclease-free centrifuge tube, avoiding the pellet and beads.

**4. Buffer Addition:**

- Add **200 µL** of **Buffer GB** to the centrifuge tube, mix thoroughly, and incubate at 70°C for 10 minutes.

**5. Ethanol Addition:**

- Add **200 µL** of **100% ethanol** to the centrifuge tube, mix well (precipitate may form after adding ethanol), and then centrifuge at 12,000 x g for 2 minutes.

**6. Supernatant Transfer:**

- Place the **DNA Spin Columns** into the **Collection Tubes** and transfer all the mixture to the DNA Spin Columns, avoiding the pellet.

**7. DNA Binding via Initial Centrifugation:**

- Centrifuge at 12,000 x g for 30 seconds.
- Discard the flow-through and return the DNA Spin Columns to the Collection Tube.

**8. Wash Steps:**

- Add **500 µL** of **Buffer PD** to the DNA Spin Columns, centrifuge at 12,000 x g for 30 seconds, discard the flow-through, and return the DNA Spin Columns to the Collection Tube.
- Add **600 µL** of **Buffer PW** to the DNA Spin Columns, centrifuge at 12,000 x g for 30 seconds, discard the flow-through, and return the DNA Spin Columns to the Collection Tube.
- **Repeat** the Buffer PW wash step.

**9. Final Centrifugation:**

- Centrifuge at 12,000 x g for 2 minutes to remove any residue.

**10. Elution:**

- Place the DNA Spin Columns into a **new 1.5 mL centrifuge tube** and let sample stand at room temperature for 5-10 minutes.
- Add **50-100 µL of Buffer TE** or nuclease-free water to the center of the membrane (note: pre-heating the Buffer TE or nuclease-free water to 65°C can improve elution efficiency).
- Centrifuge at 12,000 x g for 2 minutes to collect DNA.
- Optional: for increased yield, re-add the flow-through to the center of the membrane, let it stand for 5 minutes at room temperature, and centrifuge again at 12,000 x g for 2 minutes.

**11. Storage:**

- Store the purified genomic at -80°C for long-term storage. Avoid repeated freezing and thawing of the DNA.

**NOTES:**

1. **DISCLAIMER:** TO THE EXTENT ALLOWED BY LAW, MEDIRES CORP. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.
2. **Important Licensing Information:** These products may be covered by one or more Limited Use Label Licenses. By use of this product, you accept the terms and conditions of all applicable limited Use Label Licenses.
3. ©2024 MEDIRES CORP. All rights reserved. All trademarks are the property of MEDIRES CORP. and its subsidiaries.
4. For Research Use Only. Not for use in diagnostic procedures.